


Interim Method for the Determination of
Asbestos in Bulk Insulation Samples

Research Triangle Inst.
Research Triangle Park, NC

Prepared for

Environmental Monitoring Systems Lab.
Research Triangle Park, NC

Mar 82

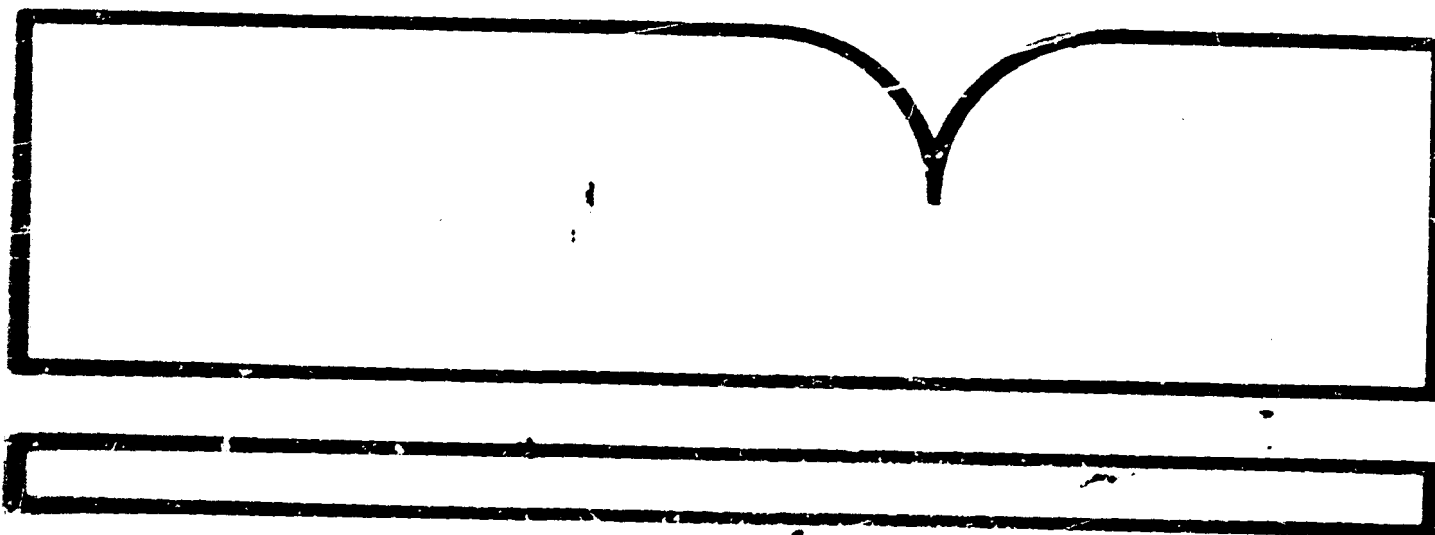

INTERIM METHOD FOR THE DETERMINATION OF ASBESTOS IN BULK INSULATION SAMPLES

by

D. E. Lentzen
E. P. Brantly, Jr.
K. W. Gold
L. E. Myers

Research Triangle Institute
P.O. Box 12194
Research Triangle Park, NC 27709

Contract No. 68-02-3431

Project Officer: J. J. Breen
Office of Pesticides and
Toxic Substances
U.S. Environmental Protection Agency
Washington, DC
20460

Project Officer: M. E. Beard
Environmental Monitoring Systems
Laboratory
U.S. Environmental Protection Agency
Research Triangle Park, NC
27711

TECHNICAL REPORT DATA		
<i>(Please read instructions on the reverse before completing)</i>		
1. REPORT NO. EPA-600/4-82-020	2. ORD Report	3. RECIPIENT'S ACCESSION NO. PB83 153643
4. TITLE AND SUBTITLE INTERIM METHOD FOR THE DETERMINATION OF ASBESTOS IN BULK INSULATION SAMPLES	5. REPORT DATE March 1982	
	6. PERFORMING ORGANIZATION CODE 43U-2069-74	
7. AUTHOR(S) D.E. Lentzen, E.P. Brantly, Jr., K.W. Gold, and L.E. Myers	8. PERFORMING ORGANIZATION REPORT NO.	
9. PERFORMING ORGANIZATION NAME AND ADDRESS Research Triangle Institute P.O. Box 12194 Research Triangle Park, N.C. 27709	10. PROGRAM ELEMENT NO.	
	11. CONTRACT GRANT NO. EPA Contract 68-02-3431	
12. SPONSORING AGENCY NAME AND ADDRESS Environmental Monitoring Systems Laboratory Research Triangle Park, NC 27711 Office of Pesticides and Toxic Substances Washington, D.C. 20460	13. TYPE OF REPORT AND PERIOD COVERED Final, January-August, 1981	
	14. SPONSORING AGENCY CODE EPA/600/06	
15. SUPPLEMENTARY NOTES		
16. ABSTRACT The U. S. Environmental Protection Agency Asbestos-in-Schools Program was established in March, 1979 to provide information and technical assistance to the public for addressing problems presented by asbestos-containing insulation materials in school buildings. Because there were no existing standard procedures for the qualitative and quantitative analysis of asbestos in bulk materials, the Office of Pesticides and Toxic Substances, Washington, D.C., and the Environmental Monitoring Systems Laboratory, Research Triangle Park, NC, jointly sponsored an effort to produce a standard analytical protocol. This report presents information on the development and characterization of the standard procedures for analysis of bulk samples with polarized light microscopy (PLM) and X-Ray diffraction (XRD), and includes the <u>Interim Method for the Determination of Asbestos in Bulk Insulation Samples</u> , (October, 1981).		
17. KEY WORDS AND DOCUMENT ANALYSIS		
a. DESCRIPTORS	b. IDENTIFIERS/OPEN ENDED TERMS	c. COSATI Field/Group
18. DISTRIBUTION STATEMENT Release to Public	19. SECURITY CLASS (This Report) Unclassified	21. NO. OF PAGES 52
	20. SECURITY CLASS (This page) Unclassified	22. PRICE


NOTICE

This document has been reviewed in accordance with U.S. Environmental Protection Agency policy and approved for publication. Mention of trade names or commercial products does not constitute endorsement or recommendation for use.

FOREWORD

Measurement and monitoring research efforts are designed to anticipate potential environmental problems, to support regulatory actions by developing an in-depth understanding of the nature and processes that impact health and the ecology, to provide innovative means of monitoring compliance with regulations, and to evaluate the effectiveness of health and environmental protection efforts through the monitoring of long-term trends. The Environmental Monitoring Systems Laboratory, Research Triangle Park, North Carolina, has responsibility for: assessment of environmental monitoring technology and systems; implementation of agency-wide quality assurance programs for air pollution measurement systems; and supplying technical support to other groups in the Agency including the Office of Air, Noise, and Radiation, the Office of Pesticides and Toxic Substances, and the Office of Enforcement.

This report describes the development of an Interim Method for the measurement of asbestos in bulk insulation materials. The method is designed to support the Asbestos-in-Schools Program of the Office of Pesticides and Toxic Substances.

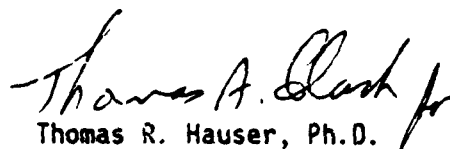
Thomas R. Hauser, Ph.D.

Director

Environmental Monitoring Systems Laboratory
Research Triangle Park, North Carolina

ABSTRACT

The U.S. Environmental Protection Agency Asbestos-in-Schools Program was established in March 1979 to provide information and technical assistance to the public for addressing problems presented by asbestos-containing insulation materials in school buildings. Because there were no existing standard procedures for the qualitative and quantitative analysis of asbestos in bulk materials, the Office of Pesticides and Toxic Substances, Washington, D.C., and the Environmental Monitoring Systems Laboratory, Research Triangle Park, NC, jointly sponsored an effort to produce a standard analytical protocol. This report presents information on the development and characterization of the standard procedures, and includes the Interim Method for the Determination of Asbestos in Bulk Insulation Samples, (October 1981).

This report is submitted in fulfillment of Contract No. 68-02-3431 by Research Triangle Institute under the sponsorship of the U.S. Environmental Protection Agency. This report covers the period January 1, 1981, to August 31, 1981, and work was completed as of October 15, 1981.

CONTENTS

	<u>Page</u>
Foreword	iii
Abstract	iv
Acknowledgment	vi
1 Introduction.	1
2 Summary	2
3 Development of the Interim Method	3
4 Overview of the Interim Method.	5
4.1 Polarized Light Microscopy.	5
4.2 X-Ray Powder Diffraction	6
References	9
Appendix	
Interim Method for the Determination of Asbestos in Bulk Insulation Samples (October 1981)	10

ACKNOWLEDGMENT

This method is based in part on contributions and review comments of participants in the symposium "Methods Definition for the Polarized Light Microscope and X-Ray Diffraction Analysis of Bulk Samples for Asbestos," U.S. Bureau of Mines, Avondale Research Center, Avondale, Maryland, October 23-24, 1979.

SECTION 1

INTRODUCTION

In March 1979, the Environmental Protection Agency (EPA) established the Asbestos-in-Schools Program to provide information and technical assistance to the public for addressing the problems presented by asbestos-containing materials in school buildings (School Asbestos Program, 44 FR 54676). Technical information concerning the identification and control of potential exposures to asbestos fibers released from friable insulation materials is available to states and local school districts. Information and assistance are available through several channels, including EPA-designated Regional Asbestos Coordinators, toll-free telephone numbers maintained by the Research Triangle Institute and the EPA Office of Pesticides and Toxic Substances (OPTS) Industry Assistance Office, and several guidance documents and program reports published by EPA.^{1 2 3}

Because there were no existing standard procedures for the analysis of asbestos in bulk materials, the Office of Pesticides and Toxic Substances, Washington, D.C., and the Environmental Monitoring Systems Laboratory, Research Triangle Park, NC, jointly sponsored an effort to produce a practical and objective analytical protocol. The Interim Method for the Determination of Asbestos in Bulk Insulation Samples has been developed by the Research Triangle Institute, on contract to EMSL and OPTS. The Interim Method includes procedures for qualitative and quantitative analysis of bulk samples by polarized light microscopy (PLM) and X-ray powder diffraction (XRD). Limited characterization of the method has been obtained in an interlaboratory study, the results of which may be found in the EPA report Bulk Sample Analysis for Asbestos Content: Evaluation of the Tentative Method (October 1981). Revisions to the original draft of the method based on the results of the evaluation testing have been incorporated into this document. Further testing is anticipated; therefore, the appended procedures are presented as an Interim Method and are subject to revision.

SECTION 2

SUMMARY

An Interim Method has been developed for the determination of asbestos in bulk samples by polarized light microscopy (PLM) and X-ray powder diffraction (XRD). EPA sponsored development of the method as part of its Asbestos-in-Schools Program.

Initial drafts of the method were based on information presented in a symposium attended by representatives of public, private, and university laboratories involved in the analysis of asbestos. Following review of the draft by symposium attendees, an interlaboratory study was conducted to collect data for the preliminary characterization of the method. Revisions to the method were made pursuant to results of the study.

The Interim Method for the Determination of Asbestos in Bulk Insulation Samples (October 1981) includes protocols for the qualitative and quantitative analysis of bulk samples by PLM and XRD, and is appended to this report.

SECTION 3

DEVELOPMENT OF THE INTERIM METHOD

The Environmental Protection Agency recommended the use of polarized light microscopy (PLM) and X-ray powder diffraction (XRD) for the analysis of bulk samples in its initial guidance document¹ for the Asbestos-in-Schools Program. PLM was recommended as the primary analytical technique, to be supplemented by XRD as necessary. Part I of the guidance document included guidelines for PLM and XRD analysis, but noted the lack of a standard protocol for bulk sample analysis. OPTS and EMSL asked the Research Triangle Institute to develop a standard protocol for the qualitative and quantitative analysis of asbestos in bulk samples.

The task was initiated by reviewing existing literature and conferring with persons knowledgeable in techniques of mineral analysis. This initial survey supported the finding that PLM and XRD were currently the most appropriate methods for use in the asbestos abatement program.

A symposium was organized for the discussion of techniques for sample preparation and qualitative and quantitative analysis by PLM and XRD. Representatives of the Federal government and analysts from public, private, and university laboratories attended the symposium "Methods Definition for the Polarized Light Microscope and X-Ray Diffraction Analysis of Bulk Samples for Asbestos" held at the U.S. Bureau of Mines, Avondale Research Center, Avondale, Maryland, October 23-24, 1979. Information presented at the symposium was used in the preparation of an initial draft of the method.

The draft protocols were distributed to symposium participants and other interested parties for review. Revisions were made to the drafts based on comments received in two rounds of review, resulting in the Tentative Method for the Determination of Asbestiform Minerals in Bulk Insulation Samples (March 1980).

An interlaboratory study was conducted for the evaluation of the Tentative Method. Results of the study are reported in the EPA report

Bulk Sample Analysis for Asbestos Content: Evaluation of the Tentative Method (October 1981). Results of the interlaboratory study and comments received from participating laboratories have been used to further revise the protocols. Because data collected in the evaluation study indicate the need for additional testing, the protocols have been designated an Interim Method and are subject to revision. The Interim Method for the Determination of Asbestos in Bulk Insulation Samples (October 1981) is appended hereto.

SECTION 4

OVERVIEW OF THE INTERIM METHOD

4.1 POLARIZED LIGHT MICROSCOPY

The PLM protocol specifies the use of classical petrographic techniques for the identification of asbestos fibers. Subsamples of bulk materials are prepared by appropriate techniques, immersed in an oil of known refractive index, and examined with a light microscope using both single and crossed polarizing filters. Asbestos fibers are positively identified by the observation of six optical properties: morphology, color and pleochroism, refractive indices (or dispersion staining colors), birefringence, extinction characteristics, and sign of elongation.

There are several deterrents to the reliable quantitation of asbestos in bulk samples by PLM, including the variability of interferences and matrices, the small amount of sample examined, and variation in the optical properties of asbestos minerals. Additionally, it is somewhat difficult to apply the results of analysis by optical microscopy to the current Federal regulations. In April 1973, EPA banned the spray application of insulation products containing more than 1 percent asbestos by weight (38 FR 2820). PLM methods measure the relative area occupied by asbestos fiber and matrix material within the microscope fields of view. The relationship between the percent asbestos of a sample on a weight basis and the percent asbestos by area as determined by PLM analysis cannot be reliably characterized for bulk samples in general.

Quantitation is performed in the PLM procedure by a point counting technique. Point sampling is used in various fields of study to determine the relative area within specified boundaries occupied by a particular subject (type of rock, soil, plant, etc.). Point counting is used in petrography to determine relative areas of minerals in thin sections of rock. The technique assumes that particles within the field of view are of

equal thickness and are randomly oriented with respect to the microscope light path.⁴ Testing has indicated that despite the violation of these assumptions, point counting may be applied to the quantitation of asbestos in bulk samples with less between-laboratory variability than the more subjective estimation techniques currently in use.

Information that allows preliminary characterization of the PLM protocol was obtained in an interlaboratory study.⁵ Samples analyzed in the study consisted of a single type of asbestos, either chrysotile or amosite, in a relatively simple matrix consisting primarily of gypsum. The results of quantitative analysis using the point counting procedure were more accurate than those produced by other methods of PLM quantitation. This was particularly true when PLM results were assumed to be a direct measure of the amount of asbestos present (i.e., ignoring the problem of relating PLM area percent results with the known weight percent values). Under the same assumption, data produced by point counting and that produced by other methods were similarly precise.

The relationship between the reported PLM data and the known values was investigated with linear regression in natural logarithmic coordinates. The principle sources of variance in the area-weight relationship were differences among laboratories and differences between asbestos types (chrysotile and amosite). When variance due to these sources was removed, the residual error of point counting data was less than that of data by other quantitation procedures.

Sources of between-laboratory variance include differences in sample preparation procedures, subsampling techniques, application of the point counting procedure, quality of optical instruments, and level of training and expertise of the analyst. Research is in progress on ways in which the application of PLM procedures may be further standardized and the accuracy of PLM data improved.

4.2 X-RAY POWDER DIFFRACTION

The XRD protocol specifies procedures for both the qualitative and quantitative analysis of asbestos in bulk materials. Since XRD affords information only on crystal lattice structure and not on gross crystal

morphology, this technique cannot distinguish between the asbestos minerals and their nonasbestiform varieties. Particle morphologies must be determined by an optical technique such as PLM. Therefore, it is recommended that the XRD method be used as a corroborative procedure with PLM and not as an independent analytical method.

Qualitative analysis of bulk materials by XRD is performed with a minimum of matrix reduction. Samples are initially scanned over limited diagnostic peak regions for the serpentine and amphibole minerals, using standard slow-scanning methods for bulk sample analysis. All samples that exhibit diffraction peaks in the diagnostic peak regions for asbestiform minerals are submitted to a full qualitative XRD scan. Sample constituents are identified by comparison of the sample diffraction pattern with standard reference powder diffraction patterns.⁶ When subsequent quantitation is required, particular note is made of possible interferences.

The thin-layer procedure for quantitation of asbestos in bulk samples by XRD is a modification of the NIOSH-recommended method for the analysis of chrysotile in air samples.⁷ This procedure involves initial comminution of the bulk material to approximately 10 μm by cryogenic milling techniques and deposition of an accurately known amount of the sample on a silver membrane filter. The mass of asbestos minerals is determined by measuring the integrated area of the selected diagnostic peak, correcting for matrix absorption effects, and comparing with suitable external standards. Although there is ample evidence that this method is capable of measuring microgram quantities of asbestos in relatively simple systems with reasonable accuracy, precision, and speed,⁸ its reliability for quantitative analysis of asbestos minerals in bulk samples has not been fully characterized.

Analytical problems and limitations of the method are clearly identified in the protocol. These include problems of sample preparation, particle size and preferred orientation effects, matrix interference and absorption effects, intraspecies interference by nonasbestos varieties of the asbestos minerals, and problems associated with assuring that reference standards are comparable to samples in chemical composition, particle size distribution, and crystallinity.

A relatively small number of laboratories provided XRD data in the interlaboratory study. None of the laboratories that participated in the

study followed the XRD protocol exactly. Some laboratories used semiquantitative thick-layer or bulk methods and others used variations of the thin-layer method. Because of the small number of laboratories that reported XRD results, and the nonequivalence of the methods employed for quantitation, it was not possible to draw any firm conclusions about the accuracy and precision of the XRD protocol. However, the results of the study did indicate that the bulk methods of analysis were at least as precise and accurate as the more rigorous thin-layer methods as applied by participating laboratories.

Recent reports in the literature have also indicated that bulk methods of analysis may provide reliable quantitative analysis of asbestos at low concentrations in a variety of matrices.^{9 10} Since XRD analysis for asbestos by bulk methods would be much less difficult and costly than thin-layer methods, further research is being directed toward comparing these methods and assessing their potential for satisfying current EPA requirements. At this time, it appears that semiquantitative bulk methods of analysis for asbestos may be acceptable as an alternative to the thin-layer method of analysis detailed in the XRD protocol.

REFERENCES

1. U.S. Environmental Protection Agency, Asbestos-Containing Materials in School Buildings: A Guidance Document, Parts 1 and 2, EPA/OTS No. C00090, March 1979.
2. D. Lucas, T. Hartwell, and A. V. Rao, Asbestos-Containing Materials in School Buildings: Guidance for Asbestos Analytical Programs, EPA 560/13-80-017A, U.S. Environmental Protection Agency, December 1980, 96 pp.
3. E. P. Brantly, Jr., and D. E. Lentzen, Asbestos-Containing Materials in School Buildings: Bulk Sample Analysis Quality Assurance Program, EPA 560/13-80-23, U.S. Environmental Protection Agency, August 1980, 27 pp.
4. F. Chayes, Petrographic Modal Analysis: An Elementary Statistical Appraisal, New York: John Wiley & Sons, 1956, 133 pp.
5. E. P. Brantly, Jr., K. W. Gold, L. E. Myers, and D. E. Lentzen, Bulk Sample Analysis for Asbestos Content: Evaluation of the Tentative Method, U.S. Environmental Protection Agency, October 1981.
6. JCPDS-International Center for Diffraction Data Powder Diffraction File, U.S. Department of Commerce, National Bureau of Standards, and Joint Committee on Powder Diffraction Studies, Swarthmore, Pennsylvania.
7. NIOSH Manual of Analytical Methods, Volume 5, U.S. Department of Health, Education and Welfare, August 1979, pp. 309-1 to 309-9.
8. B. A. Lange and J. C. Haartz, Determination of microgram quantities of asbestos by X-ray diffraction: Chrysotile in thin dust layers of matrix material, Anal. Chem., 51(4):520-525, 1979.
9. H. Dunn and J. H. Stewart, Jr., Quantitative determination of chrysotile in building materials, The Microscope, 29(1):39-45, 1981.
10. M. Taylor, Methods for quantitative determination of asbestos and quartz in bulk samples using X-ray diffraction, The Analyst, 103(1231):1009-1020, 1978.

APPENDIX
INTERIM METHOD FOR THE DETERMINATION OF ASBESTOS
IN BULK INSULATION SAMPLES

INTERIM METHOD FOR THE DETERMINATION OF ASBESTOS
IN BULK INSULATION SAMPLES

An interim method is carefully drafted from available source information. This method is still under investigation and therefore is subject to revision.

Prepared for

Office of Pesticides and
Toxic Substances
U.S. Environmental Protection
Agency
Washington, DC
20460

Environmental Monitoring
Systems Laboratory
U.S. Environmental Protection
Agency
Research Triangle Park, NC
27711

October 1981

SECTION 1

POLARIZED LIGHT MICROSCOPY

1.1 PRINCIPLE AND APPLICABILITY

Bulk samples of building materials taken for asbestos identification are first examined for homogeneity and preliminary fiber identification at low magnification. Positive identification of suspect fibers is made by analysis of subsamples with the polarized light microscope.

The principles of optical mineralogy are well established.^{1 2} A light microscope equipped with two polarizing filters is used to observe specific optical characteristics of a sample. The use of plane polarized light allows the determination of refractive indices along specific crystallographic axes. Morphology and color are also observed. A retardation plate is placed in the polarized light path for determination of the sign of elongation using orthoscopic illumination. Orientation of the two filters such that their vibration planes are perpendicular (crossed polars) allows observation of the birefringence and extinction characteristics of anisotropic particles.

Quantitative analysis involves the use of point counting. Point counting is a standard technique in petrography for determining the relative areas occupied by separate minerals in thin sections of rock. Background information on the use of point counting² and the interpretation of point count data³ is available.

This method is applicable to all bulk samples of friable insulation materials submitted for identification and quantitation of asbestos components.

1.2 RANGE

The point counting method may be used for analysis of samples containing from 0 to 100 percent asbestos. The upper detection limit is 100 percent. The lower detection limit is less than 1 percent.

1.3 INTERFERENCES

Fibrous organic and inorganic constituents of bulk samples may interfere with the identification and quantitation of the asbestos mineral content. Spray-on binder materials may coat fibers and affect color or obscure optical characteristics to the extent of masking fiber identity. Fine particles of other materials may also adhere to fibers to an extent sufficient to cause confusion in identification. Procedures that may be used for the removal of interferences are presented in Section 1.7.2.2.

1.4 PRECISION AND ACCURACY

Adequate data for measuring the accuracy and precision of the method for samples with various matrices are not currently available. Data obtained for samples containing a single asbestos type in a simple matrix are available in the EPA report Bulk Sample Analysis for Asbestos Content: Evaluation of the Tentative Method.⁴

1.5 APPARATUS

1.5.1 Sample Analysis

A low-power binocular microscope, preferably stereoscopic, is used to examine the bulk insulation sample as received.

- Microscope: binocular, 10-45X (approximate).
- Light Source: incandescent or fluorescent.
- Forceps, Dissecting Needles, and Probes
- Glassine Paper or Clean Glass Plate

Compound microscope requirements: A polarized light microscope complete with polarizer, analyzer, port for wave retardation plate, 360° graduated rotating stage, substage condenser, lamp, and lamp iris.

- Polarized Light Microscope: described above.
- Objective Lenses: 10X, 20X, and 40X or near equivalent.
- Dispersion Staining Objective Lens (optional)
- Ocular Lens: 10X minimum.
- Eyepiece Reticle: cross hair or 25 point Chalkley Point Array.
- Compensator Plate: 550 millimicron retardation.

1.5.2 Sample Preparation

Sample preparation apparatus requirements will depend upon the type of insulation sample under consideration. Various physical and/or chemical means may be employed for an adequate sample assessment.

- Ventilated Hood or negative pressure glove box.
- Microscope Slides
- Coverslips
- Mortar and Pestle: agate or porcelain. (optional)
- Wylie Mill (optional)
- Beakers and Assorted Glassware (optional)
- Centrifuge (optional)
- Filtration apparatus (optional)
- Low temperature asher (optional)

1.6 REAGENTS

1.6.1 Sample Preparation

- Distilled Water (optional)
- Dilute CH₃COOH: ACS reagent grade (optional)
- Dilute HCl: ACS reagent grade (optional)
- Sodium metaphosphate (NaPO₃)₆ (optional)

1.6.2 Analytical Reagents

- Refractive Index Liquids: 1.490-1.570, 1.590-1.720 in increments of 0.002 or 0.004.
- Refractive Index Liquids for Dispersion Staining: high-dispersion series, 1.550, 1.605, 1.630 (optional).
- UICC Asbestos Reference Sample Set: Available from: UICC MRC Pneumoconiosis Unit, Llandough Hospital, Penarth, Glamorgan CF6 1XW, UK, and commercial distributors.
- Tremolite-asbestos (source to be determined)
- Actinolite-asbestos (source to be determined)

1.7 PROCEDURES

NOTE: EXPOSURE TO AIRBORNE ASBESTOS FIBERS IS A HEALTH HAZARD. BULK SAMPLES SUBMITTED FOR ANALYSIS ARE USUALLY FRIABLE AND MAY RELEASE FIBERS DURING HANDLING OR MATRIX REDUCTION STEPS. ALL SAMPLE AND SLIDE PREPARATIONS

SHOULD BE CARRIED OUT IN A VENTILATED HOOD OR GLOVE BOX WITH CONTINUOUS AIRFLOW (NEGATIVE PRESSURE). HANDLING OF SAMPLES WITHOUT THESE PRECAUTIONS MAY RESULT IN EXPOSURE OF THE ANALYST AND CONTAMINATION OF SAMPLES BY AIRBORNE FIBERS.

1.7.1 Sampling

Samples for analysis of asbestos content shall be taken in the manner prescribed in Reference 5 and information on design of sampling and analysis programs may be found in Reference 6. If there are any questions about the representative nature of the sample, another sample should be requested before proceeding with the analysis.

1.7.2 Analysis

1.7.2.1 Gross Examination--

Bulk samples of building materials taken for the identification and quantitation of asbestos are first examined for homogeneity at low magnification with the aid of a stereomicroscope. The core sample may be examined in its container or carefully removed from the container onto a glassine transfer paper or clean glass plate. If possible, note is made of the top and bottom orientation. When discrete strata are identified, each is treated as a separate material so that fibers are first identified and quantified in that layer only, and then the results for each layer are combined to yield an estimate of asbestos content for the whole sample.

1.7.2.2 Sample Preparation--

Bulk materials submitted for asbestos analysis involve a wide variety of matrix materials. Representative subsamples may not be readily obtainable by simple means in heterogeneous materials, and various steps may be required to alleviate the difficulties encountered. In most cases, however, the best preparation is made by using forceps to sample at several places from the bulk material. Forcep samples are immersed in a refractive index liquid on a microscope slide, teased apart, covered with a cover glass, and observed with the polarized light microscope.

Alternatively, attempts may be made to homogenize the sample or eliminate interferences before further characterization. The selection of appropriate procedures is dependent upon the samples encountered and personal preference. The following are presented as possible sample preparation steps.

A mortar and pestle can sometimes be used in the size reduction of soft or loosely bound materials though this may cause matting of some samples. Such samples may be reduced in a Wylie mill. Apparatus should be clean and extreme care exercised to avoid cross-contamination of samples. Periodic checks of the particle sizes should be made during the grinding operation so as to preserve any fiber bundles present in an identifiable form. These procedures are not recommended for samples that contain amphibole minerals or vermiculite. Grinding of amphiboles may result in the separation of fiber bundles or the production of cleavage fragments with aspect ratios greater than 3:1. Grinding of vermiculite may also produce fragments with aspect ratios greater than 3:1.

Acid treatment may occasionally be required to eliminate interferences. Calcium carbonate, gypsum, and bassanite (plaster) are frequently present in sprayed or trowelled insulations. These materials may be removed by treatment with warm dilute acetic acid. Warm dilute hydrochloric acid may also be used to remove the above materials. If acid treatment is required, wash the sample at least twice with distilled water, being careful not to lose the particulates during decanting steps. Centrifugation or filtration of the suspension will prevent significant fiber loss. The pore size of the filter should be 0.45 micron or less. Caution: prolonged acid contact with the sample may alter the optical characteristics of chrysotile fibers and should be avoided.

Coatings and binding materials adhering to fiber surfaces may also be removed by treatment with sodium metaphosphate.⁷ Add 10 mL of 10 g/L sodium metaphosphate solution to a small (0.1 to 0.5 mL) sample of bulk material in a 15-mL glass centrifuge tube. For approximately 15 seconds each, stir the mixture on a vortex mixer, place in an ultrasonic bath and then shake by hand. Repeat the series. Collect the dispersed solids by centrifugation at 1000 rpm for 5 minutes. Wash the sample three times by suspending in 10 mL distilled water and recentrifuging. After washing, resuspend the pellet in 5 mL distilled water, place a drop of the suspension on a microscope slide, and dry the slide at 110° C.

In samples with a large portion of cellulosic or other organic fibers, it may be useful to ash part of the sample and view the residue. Ashing should be performed in a low temperature asher. Ashing may also be performed in a muffle furnace at temperatures of 500° C or lower. Temperatures of 550° C or higher will cause dehydroxylation of the asbestos minerals, resulting in changes

of the refractive index and other key parameters. If a muffle furnace is to be used, the furnace thermostat should be checked and calibrated to ensure that samples will not be heated at temperatures greater than 550° C.

Ashing and acid treatment of samples should not be used as standard procedures. In order to monitor possible changes in fiber characteristics, the material should be viewed microscopically before and after any sample preparation procedure. Use of these procedures on samples to be used for quantitation requires a correction for percent weight loss.

1.7.2.3 Fiber Identification--

Positive identification of asbestos requires the determination of the following optical properties.

- Morphology
- Color and pleochroism
- Refractive indices
- Birefringence
- Extinction characteristics
- Sign of elongation

Table 1-1 lists the above properties for commercial asbestos fibers. Figure 1-1 presents a flow diagram of the examination procedure. Natural variations in the conditions under which deposits of asbestiform minerals are formed will occasionally produce exceptions to the published values and differences from the UICC standards. The sign of elongation is determined by use of the compensator plate and crossed polars. Refractive indices may be determined by the Becke line test. Alternatively, dispersion staining may be used. Inexperienced operators may find that the dispersion staining technique is more easily learned, and should consult Reference 9 for guidance. Central stop dispersion staining colors are presented in Table 1-2. Available high-dispersion (HD) liquids should be used.

1.7.2.4 Quantitation of Asbestos Content--

Asbestos quantitation is performed by a point-counting procedure. An ocular reticle (cross-hair or point array) is used to visually superimpose a point or points on the microscope field of view. Record the number of points positioned directly above each kind of particle or fiber of interest. Score only points directly over asbestos fibers or nonasbestos matrix material. Do

TABLE 1-1. OPTICAL PROPERTIES OF ASBESTOS FIBERS

Mineral	Morphology, color ^a	Refractive indices ^b		Birefringence	Extinction	Sign of elongation
		α	γ			
Chrysotile (asbestiform serpentine)	Wavy fibers. Fiber bundles have splayed ends and "kinks". Aspect ratio typically >10:1. Colorless, nonpleochroic.	1.493-1.560	1.517-1.562 ^f (normally 1.556)	.008	to fiber length	+ (length slow)
Amosite (asbestiform grunerite)	Straight, rigid fibers. Aspect ratio typically >10:1. Colorless to brown, nonpleochroic or weakly so. Opaque inclusions may be present.	1.635-1.696	1.655-1.729 ^f (normally 1.696-1.710)	.020-.033	to fiber length	+ (length slow)
Crocidolite (asbestiform riebeckite)	Straight, rigid fibers. Thick fibers and bundles common, blue to purple-blue in color. Pleochroic. Birefringence is generally masked by blue color.	1.654-1.701	1.668-1.717 ^e (normally close to 1.700)	.014-.016	to fiber length	- (length: fast)
Anthophyllite- asbestos	Straight fibers and acicular cleavage fragments. Some composite fibers. Aspect ratio <10:1. Colorless to light brown.	1.596-1.652	1.615-1.676 ^f	.019-.024	to fiber length	+ (length slow)
Tremolite- actinolite- asbestos	Normally present as acicular or prismatic cleavage fragments. Single crystals predominate, aspect ratio <10:1. Colorless to pale green.	1.599-1.668	1.622-1.688 ^f	.023-.020	Oblique extinction, 10-20° for fragments. Composite fibers show extinction.	+ (length slow)

^aFrom reference 5; colors cited are seen by observation with plane polarized light.

^bFrom references 5 and 8.

^cFibers subjected to heating may be brownish.

^dFibers defined as having aspect ratio >3:1.

^e|| to fiber length.

^f|| to fiber length.

Polarized light microscopy analysis: For each type of material identified by examination of sample at low magnification. Mount spacially dispersed sample in 1.550 RI liquid. (If using dispersion staining, mount in 1.550 HD.) View at 100X with both plane polarized light and crossed polars. More than one fiber type may be present.

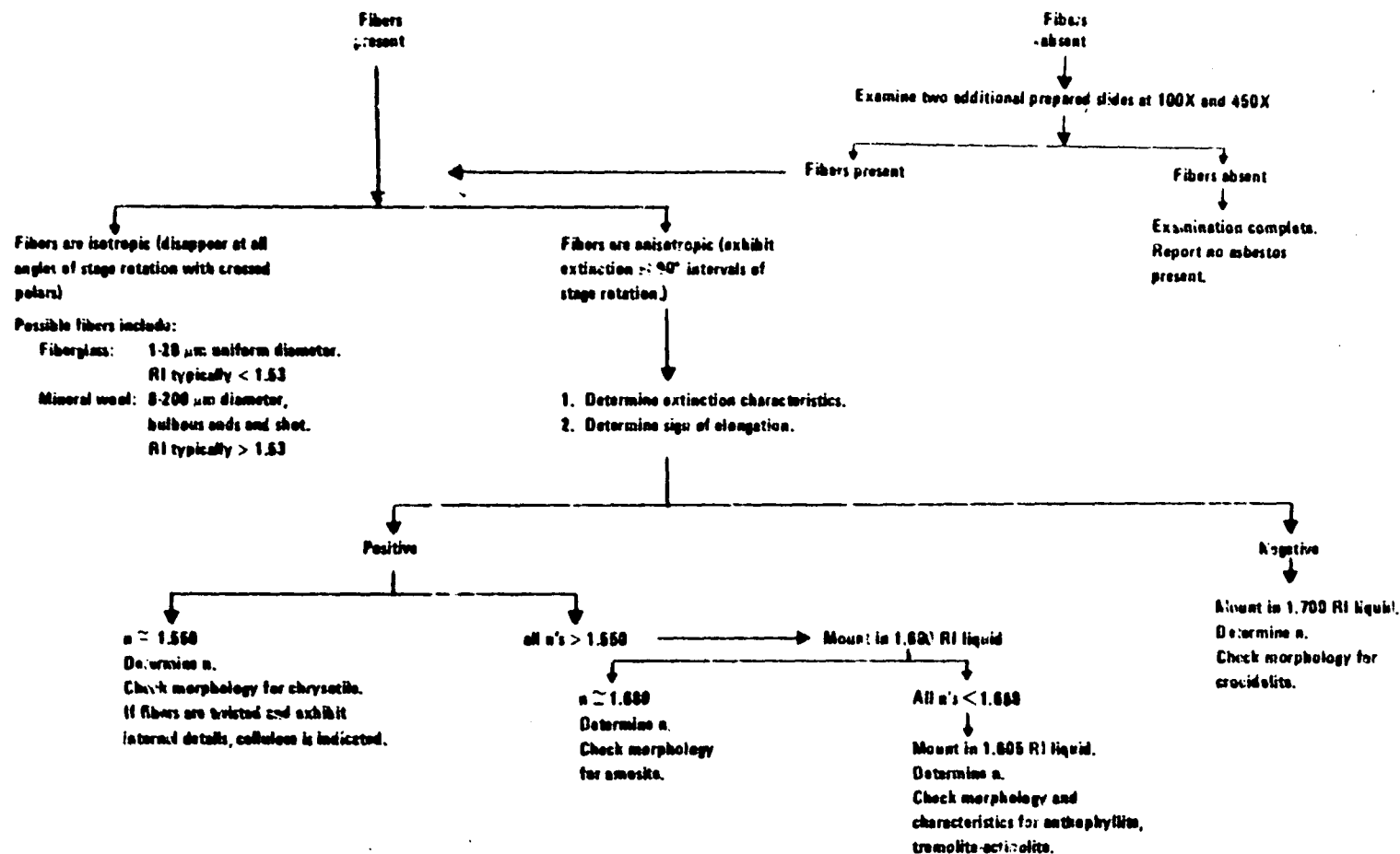


Figure 1-1. Flow chart for analysis of bulk samples by polarized light microscopy.

TABLE 1-2. CENTRAL STOP DISPERSION STAINING COLORS^a

Mineral	RI Liquid	n_{\perp}	n_{\parallel}
Chrysotile	1.550 ^{HD}	Blue	Blue-magenta
Amosite	1.680	Blue-magenta to pale blue	Golden-yellow
	1.550 ^{HD}	Yellow to white	Yellow to white
Crocidolite ^b	1.700	Red magenta	Blue-magenta
	1.550 ^{HD}	Yellow to white	Yellow to white
Anthophyllite	1.605 ^{HD}	Blue	Gold to gold-magenta
Tremolite	1.605 ^{HD^c}	Pale blue	Yellow
Actinolite	1.605 ^{HD}	Gold-magenta to blue	Gold
	1.630 ^{HD^c}	Magenta	Golden-yellow

^aFrom reference 9.^bBlue absorption color.^cOblique extinction view.

not score empty points for the closest particle. If an asbestos fiber and a matrix particle overlap so that a point is superimposed on their visual intersection, a point is scored for both categories. Point counting provides a determination of the area percent asbestos. Reliable conversion of area percent to percent of dry weight is not currently feasible unless the specific gravities and relative volumes of the materials are known.

For the purpose of this method, "asbestos fibers" are defined as having an aspect ratio greater than 3:1 and being positively identified as one of the minerals in Table 1-1.

A total of 400 points superimposed on either asbestos fibers or nonasbestos matrix material must be counted over at least eight different preparations of representative subsamples. Take eight forcep samples and mount each separately with the appropriate refractive index liquid. The preparation should not be heavily loaded. The sample should be uniformly dispersed to avoid overlapping particles and allow 25-50 percent empty area within the fields of view. Count 50 nonempty points on each preparation, using either

- A cross-hair reticle and mechanical stage; or
- A reticle with 25 points (Chalkley Point Array) and counting at least 2 randomly selected fields.

For samples with mixtures of isotropic and anisotropic materials present, viewing the sample with slightly uncrossed polars or the addition of the compensator plate to the polarized light path will allow simultaneous discrimination of both particle types. Quantitation should be performed at 100X or at the lowest magnification of the polarized light microscope that can effectively distinguish the sample components. Confirmation of the quantitation result by a second analyst on some percentage of analyzed samples should be used as standard quality control procedure.

The percent asbestos is calculated as follows:

$$\% \text{ asbestos} = (a/n) 100\%$$

where

a = number of asbestos counts,

n = number of nonempty points counted (400).

If $a = 0$, report "No asbestos detected." If $0 < a \leq 3$, report "<1% asbestos".

The value reported should be rounded to the nearest percent.

1.8 REFERENCES

1. Paul F. Kerr, Optical Mineralogy, 4th ed., New York, McGraw-Hill, 1977.
2. E. M. Chamot and C. W. Mason, Handbook of Chemical Microscopy, Volume One, 3rd ed., New York: John Wiley & Sons, 1958.
3. F. Chayes, Petrographic Modal Analysis: An Elementary Statistical Appraisal, New York: John Wiley & Sons, 1956.
4. E. P. Brantly, Jr., K. W. Gold, L. E. Myers, and D. E. Lentzen, Bulk Sample Analysis for Asbestos Content: Evaluation of the Tentative Method, U.S. Environmental Protection Agency, October 1981.
5. U.S. Environmental Protection Agency, Asbestos-Containing Materials in School Buildings: A Guidance Document, Parts 1 and 2, EPA/OTS No. C00090, March 1979.
6. D. Lucas, T. Hartwell, and A. V. Rao, Asbestos-Containing Materials in School Buildings: Guidance for Asbestos Analytical Programs, EPA 560/13-80-017A, U.S. Environmental Protection Agency, December 1980, 96 pp.
7. D. H. Taylor and J. S. Bloom, Hexametaphosphate pretreatment of insulation samples for identification of fibrous constituents, Microscope, 28, 1980.
8. W. J. Campbell, R. L. Blake, L. L. Brown, E. E. Cather, and J. J. Sjöberg, Selected Silicate Minerals and Their Asbestiform Varieties: Mineralogical Definitions and Identification-Characterization, U.S. Bureau of Mines Information Circular 8751, 1977.
9. Walter C. McCrone, Asbestos Particle Atlas, Ann Arbor: Ann Arbor Science Publishers, June 1980.

SECTION 2

X-RAY POWDER DIFFRACTION

2.1 PRINCIPLE AND APPLICABILITY

The principle of X-ray powder diffraction (XRD) analysis is well established.^{1 2} Any solid, crystalline material will diffract an impinging beam of parallel, monochromatic X-rays whenever Bragg's Law,

$$\lambda = 2d \sin \theta,$$

is satisfied for a particular set of planes in the crystal lattice, where

λ = the X-ray wavelength, Å;

d = the interplanar spacings of the set of reflecting lattice planes, Å; and

θ = the angle of incidence between the X-ray beam and the reflecting lattice planes.

By appropriate orientation of a sample relative to the incident X-ray beam, a diffraction pattern can be generated that, in most cases, will be uniquely characteristic of both the chemical composition and structure of the crystalline phases present.

Unlike optical methods of analysis, however, XRD cannot determine crystal morphology. Therefore, in asbestos analysis, XRD does not distinguish between fibrous and nonfibrous forms of the serpentine and amphibole minerals (Table 2-1). However, when used in conjunction with optical methods such as polarized light microscopy (PLM), XRD techniques can provide a reliable analytical method for the identification and characterization of asbestiform minerals in bulk materials.

For qualitative analysis by XRD methods, samples are initially scanned over limited diagnostic peak regions for the serpentine (~7.4 Å) and amphibole (8.2-8.5 Å) minerals (Table 2-2). Standard slow-scanning methods for bulk sample analysis may be used for materials shown by PLM to contain significant amounts of asbestos (>5-10 percent). Detection of minor or trace amounts of

TABLE 2-1. THE ASBESTOS MINERALS AND THEIR
NONASBESTIFORM ANALOGS

Asbestiform	Nonasbestiform
SERPENTINE	
Chrysotile	Antigorite, lizardite
AMPHIBOLE	
Anthophyllite asbestos	Anthophyllite
Cummingtonite-grunerite asbestos ("Amosite")	Cummingtonite-grunerite
Crocidolite	Riebeckite
Tremolite asbestos	Tremolite
Actinolite asbestos	Actinolite

TABLE 2-2. PRINCIPAL LATTICE SPACINGS OF ASBESTIFORM MINERALS^a

Minerals	Principal d-spacings (Å) and relative intensities			JCPDS Powder diffraction file ³ number
Chrysotile	7.37 ₁₀₀	3.55 ₇₀	4.57 ₅₀	21-543 ^b
	7.36 ₁₀₀	3.66 ₈₀	2.45 ₆₅	25-645
	7.10 ₁₀₀	2.33 ₈₀	3.55 ₇₀	22-1162 (theoretical)
"Amosite"	8.33 ₁₀₀	3.06 ₇₀	2.75 ₆₇₀	17-745 (nonfibrous)
	8.22 ₁₀₀	3.06 ₈₅	3.25 ₇₀	27-1170 (UICC)
Anthophyllite	3.05 ₁₀₀	3.24 ₅₀	8.26 ₅₅	9-455
	3.06 ₁₀₀	8.33 ₇₀	3.23 ₅₀	16-401 (synthetic)
Actinolite	2.72 ₁₀₀	2.54 ₁₀₀	3.40 ₈₀	25-157
Crocidolite	8.35 ₁₀₀	3.10 ₅₅	2.72 ₃₅	27-1415 (UICC)
Tremolite	8.39 ₁₀₀	3.12 ₁₀₀	2.705 ₉₀	13-437 ^b
	2.706 ₁₀₀	3.14 ₉₅	8.43 ₄₀	20-1310 ^b (synthetic)
	3.13 ₁₀₀	2.706 ₆₀	8.44 ₄₀	23-666 (synthetic mixture with richterite)

^aThis information is intended as a guide only. Complete powder diffraction data, including mineral type and source, should be referred to, to ensure comparability of sample and reference materials where possible. Additional precision XRD data on amosite, crocidolite, tremolite, and chrysotile are available from the U.S. Bureau of Mines.⁴

^bFibrosity questionable.

asbestos may require special sample preparation and step-scanning analysis. All samples that exhibit diffraction peaks in the diagnostic regions for asbestiform minerals are submitted to a full (5° - 60° 2θ ; 1° $2\theta/\text{min}$) qualitative XRD scan, and their diffraction patterns are compared with standard reference powder diffraction patterns³ to verify initial peak assignments and to identify possible matrix interferences when subsequent quantitative analysis will be performed.

Accurate quantitative analysis of asbestos in bulk samples by XRD is critically dependent on particle size distribution, crystallite size, preferred orientation and matrix absorption effects, and comparability of standard reference and sample materials. The most intense diffraction peak that has been shown to be free from interference by prior qualitative XRD analysis is selected for quantitation of each asbestiform mineral. A "thin-layer" method of analysis^{5, 6} is recommended in which, subsequent to comminution of the bulk material to $\sim 10\ \mu\text{m}$ by suitable cryogenic milling techniques, an accurately known amount of the sample is deposited on a silver membrane filter. The mass of asbestiform material is determined by measuring the integrated area of the selected diffraction peak using a step-scanning mode, correcting for matrix absorption effects, and comparing with suitable calibration standards. Alternative "thick-layer" or bulk methods,^{7, 8} may be used for semiquantitative analysis.

This XRD method is applicable as a confirmatory method for identification and quantitation of asbestos in bulk material samples that have undergone prior analysis by PLM or other optical methods.

2.2 RANGE AND SENSITIVITY

The range of the method has not been determined.

The sensitivity of the method has not been determined. It will be variable and dependent upon many factors, including matrix effects (absorption and interferences), diagnostic reflections selected, and their relative intensities.

2.3 LIMITATIONS

2.3.1 Interferences

Since the fibrous and nonfibrous forms of the serpentine and amphibole minerals (Table 2-1) are indistinguishable by XRD techniques unless special

sample preparation techniques and instrumentation are used,⁹ the presence of nonasbestiform serpentines and amphiboles in a sample will pose severe interference problems in the identification and quantitative analysis of their asbestiform analogs.

The use of XRD for identification and quantitation of asbestiform minerals in bulk samples may also be limited by the presence of other interfering materials in the sample. For naturally occurring materials the commonly associated asbestos-related mineral interferences can usually be anticipated. However, for fabricated materials the nature of the interferences may vary greatly (Table 2-3) and present more serious problems in identification and quantitation.¹⁰ Potential interferences are summarized in Table 2-4 and include the following:

- Chlorite has major peaks at 7.19 Å and 3.58 Å that interfere with both the primary (7.36 Å) and secondary (3.66 Å) peaks for chrysotile. Resolution of the primary peak to give good quantitative results may be possible when a step-scanning mode of operation is employed.
- Halloysite has a peak at 3.63 Å that interferes with the secondary (3.66 Å) peak for chrysotile.
- Kaolinite has a major peak at 7.15 Å that may interfere with the primary peak of chrysotile at 7.36 Å when present at concentrations of >10 percent. However, the secondary chrysotile peak at 3.66 Å may be used for quantitation.
- Gypsum has a major peak at 7.5 Å that overlaps the 7.36 Å peak of chrysotile when present as a major sample constituent. This may be removed by careful washing with distilled water, or by heating to 300° C to convert gypsum to plaster of paris.
- Cellulose has a broad peak that partially overlaps the secondary (3.65 Å) chrysotile peak.⁸
- Overlap of major diagnostic peaks of the amphibole asbestos minerals, amosite, anthophyllite, crocidolite, and tremolite, at approximately 8.3 Å and 3.1 Å causes mutual interference when these minerals occur in the presence of one another. In some instances, adequate resolution may be attained by using step-scanning methods and/or by decreasing the collimator slit width at the X-ray port.

TABLE 2-3. COMMON CONSTITUENTS IN INSULATION AND
WALL MATERIALS¹⁰

A. Insulation materials

Chrysotile
 "Amosite"
 Crocidolite
 *Rock wool
 *Slag wool
 *Fiber glass
 Gypsum ($\text{CaSO}_4 \cdot 2\text{H}_2\text{O}$)
 Vermiculite (micas)
 *Perlite
 Clays (kaolin)
 *Wood pulp
 *Paper fibers (talc, clay, carbonate fillers)
 Calcium silicates (synthetic)
 Opaques (chromite, magnetite inclusions in serpentine)
 Hematite (inclusions in "amosite")
 Magnesite
 *Diatomaceous earth

B. Spray finishes or paints

Bassanite
 Carbonate minerals (calcite, dolomite, vaterite)
 Talc
 Tremolite
 Anthophyllite
 Serpentine (including chrysotile)
 Amosite
 Crocidolite
 *Mineral wool
 *Rock wool
 *Slag wool
 *Fiber glass
 Clays (kaolin)
 Micas
 Chlorite
 Gypsum ($\text{CaSO}_4 \cdot 2\text{H}_2\text{O}$)
 Quartz
 *Organic binders and thickeners
 Hydromagnesite
 Wollastonite
 Opaques (chromite, magnetite inclusions in serpentine)
 Hematite (inclusions in "amosite")

*Amorphous materials--contribute only to overall scattered radiation and increased background radiation.

TABLE 2-4. INTERFERENCES IN XRD ANALYSIS
OF ASBESTIFORM MINERALS

Asbestiform mineral	Primary diagnostic peaks (approximate d-spacings, in Å)	Interference
Serpentine		
Chrysotile	7.4	Nonasbestiform serpentines (antigorite, lizardite) Chlorite Kaolinite Gypsum
	3.7	Chlorite Halloysite Cellulose
Amphibole		
"Amosite"	3.1	Nonasbestiform amphiboles (cummingtonite-grunerite, anthophyllite, riebeckite, tremolite) Mutual interferences Carbonates Talc
Anthophyllite		
Crocidolite		
Tremolite		
	8.3	Mutual interferences

- Carbonates may also interfere with quantitative analysis of the amphibole asbestos minerals, amosite, anthophyllite, crocidolite, and tremolite. Calcium carbonate (CaCO_3) has a peak at 3.035 Å that overlaps major amphibole peaks at approximately 3.1 Å when present in concentrations of >5 percent. Removal of carbonates with a dilute acid wash is possible; however, if present, chrysotile may be partially dissolved by this treatment.¹¹
- A major talc peak at 3.12 Å interferes with the primary tremolite peak at this same position and with secondary peaks of crocidolite (3.10 Å), amosite (3.06 Å), and anthophyllite (3.05 Å). In the presence of talc, the major diagnostic peak at approximately 8.3 Å should be used for quantitation of these asbestiform minerals.

The problem of intraspecies and matrix interferences is further aggravated by the variability of the silicate mineral powder diffraction patterns themselves, which often makes definitive identification of the asbestos minerals by comparison with standard reference diffraction patterns difficult. This variability results from alterations in the crystal lattice associated with differences in isomorphous substitution and degree of crystallinity. This is especially true for the amphiboles. These minerals exhibit a wide variety of very similar chemical compositions, with the result being that their diffraction patterns are characterized by having major (110) reflections of the monoclinic amphiboles and (210) reflections of the orthorhombic anthophyllite separated by less than 0.2 Å.¹²

2.3.2 Matrix Effects

If a copper X-ray source is used, the presence of iron at high concentrations in a sample will result in significant X-ray fluorescence, leading to loss of peak intensity along with increased background intensity and an overall decrease in sensitivity. This situation may be corrected by choosing an X-ray source other than copper; however, this is often accompanied both by loss of intensity and by decreased resolution of closely spaced reflections. Alternatively, use of a diffracted beam monochromator will reduce background fluorescent radiation, enabling weaker diffraction peaks to be detected.

X-ray absorption by the sample matrix will result in overall attenuation of the diffracted beam and may seriously interfere with quantitative analysis. Absorption effects may be minimized by using sufficiently "thin" samples for analysis.^{5 13 14} However, unless absorption effects are known to be the same for both samples and standards, appropriate corrections should be made by referencing diagnostic peak areas to an internal standard^{7 8} or filter substrate (Ag) peak.^{5 6}

2.3.3 Particle Size Dependence

Because the intensity of diffracted X-radiation is particle-size dependent, it is essential for accurate quantitative analysis that both sample and standard reference materials have similar particle size distributions. The optimum particle size range for quantitative analysis of asbestos by XRD has been reported to be 1 to 10 μm .¹⁵ Comparability of sample and standard reference material particle size distributions should be verified by optical microscopy (or another suitable method) prior to analysis.

2.3.4 Preferred Orientation Effects

Preferred orientation of asbestiform minerals during sample preparation often poses a serious problem in quantitative analysis by XRD. A number of techniques have been developed for reducing preferred orientation effects in "thick layer" samples.^{7 8 15} However, for "thin" samples on membrane filters, the preferred orientation effects seem to be both reproducible and favorable to enhancement of the principal diagnostic reflections of asbestos minerals, actually increasing the overall sensitivity of the method.^{12 14} (Further investigation into preferred orientation effects in both thin layer and bulk samples is required.)

2.3.5 Lack of Suitably Characterized Standard Materials

The problem of obtaining and characterizing suitable reference materials for asbestos analysis is clearly recognized. NIOSH has recently directed a major research effort toward the preparation and characterization of analytical reference materials, including asbestos standards;^{16 17} however, these are not available in large quantities for routine analysis.

In addition, the problem of ensuring the comparability of standard reference and sample materials, particularly regarding crystallite size, particle size distribution, and degree of crystallinity, has yet to be adequately addressed. For example, Langer et al.¹⁸ have observed that in insulating matrices, chrysotile tends to break open into bundles more frequently than amphiboles. This results in a line-broadening effect with a resultant decrease in sensitivity. Unless this effect is the same for both standard and sample materials, the amount of chrysotile in the sample will be underestimated by XRD analysis. To minimize this problem, it is recommended that standardized matrix reduction procedures be used for both sample and standard materials.

2.4 PRECISION AND ACCURACY

Precision of the method has not been determined.

Accuracy of the method has not been determined.

2.5 APPARATUS

2.5.1 Sample Preparation

Sample preparation apparatus requirements will depend upon the sample type under consideration and the kind of XRD analysis to be performed.

- Mortar and Pestle: Agate or porcelain.
- Razor Blades
- Sample Mill: SPEX, Inc., freezer mill or equivalent.
- Bulk Sample Holders
- Silver Membrane Filters: 25-mm diameter, 0.45- μ m pore size. Selas Corp. of America, Flotronics Div., 1957 Pioneer Road, Huntingdon Valley, PA 19006.
- Microscope Slides
- Vacuum Filtration Apparatus: Gelman No. 1107 or equivalent, and side-arm vacuum flask.
- Microbalance
- Ultrasonic Bath or Probe: Model W140, Ultrasonics, Inc., operated at a power density of approximately 0.1 W/mL, or equivalent.

- Volumetric Flasks: 1-L volume.
- Assorted Pipettes
- Pipette Bulb
- Nonserrated Forceps
- Polyethylene Wash Bottle
- Pyrex Beakers: 50-mL volume.
- Desiccator
- Filter Storage Cassettes
- Magnetic Stirring Plate and Bars
- Porcelain Crucibles
- Muffle Furnace or Low Temperature Asher

2.5.2 Sample Analysis

Sample analysis requirements include an X-ray diffraction unit, equipped with:

- Constant Potential Generator; Voltage and mA Stabilizers
- Automated Diffractometer with Step-Scanning Mode
- Copper Target X-Ray Tube: High intensity, fine focus, preferably.
- X-Ray Pulse Height Selector
- X-Ray Detector (with high voltage power supply): Scintillation or proportional counter.
- Focusing Graphite Crystal Monochromator; or Nickel Filter (if copper source is used, and iron fluorescence is not a serious problem).
- Data Output Accessories:
 - Strip Chart Recorder
 - Decade Scaler/Timer
 - Digital Printer

- Sample Spinner (optional).
- Instrument Calibration Reference Specimen: α -quartz reference crystal (Arkansas quartz standard, #180-147-00, Philips Electronics Instruments, Inc., 85 McKee Drive, Mahwah, NJ 07430) or equivalent.

2.6 REAGENTS

2.6.1 Standard Reference Materials

The reference materials listed below are intended to serve as a guide. Every attempt should be made to acquire pure reference materials that are comparable to sample materials being analyzed.

- Chrysotile: UICC Canadian, or NIEHS Plastibest. (UICC reference materials available from: UICC, MRC Pneumoconiosis Unit, Llandough Hospital, Penarth, Glamorgan, CF61XW, UK).
- Crocidolite: UICC
- Amosite: UICC
- Anthophyllite: UICC
- Tremolite Asbestos: Wards Natural Science Establishment, Rochester, N.Y.; Cyprus Research Standard, Cyprus Research, 2435 Military Ave., Los Angeles, CA 90064 (washed with dilute HCl to remove small amount of calcite impurity); India tremolite, Rajasthan State, India.
- Actinolite Asbestos

2.6.2 Adhesive

Tape, petroleum jelly, etc. (for attaching silver membrane filters to sample holders).

2.6.3 Surfactant

1 percent aerosol OT aqueous solution or equivalent.

2.6.4 Isopropanol

ACS Reagent Grade.

2.7 PROCEDURE

2.7.1 Sampling

Samples for analysis of asbestos content shall be collected as specified

in EPA Guidance Document #C0090, Asbestos-Containing Materials in School Buildings.¹⁰

2.7.2 Analysis

All samples must be analyzed initially for asbestos content by PLM. XRD should be used as an auxiliary method when a second, independent analysis is requested.

NOTE: ASBESTOS IS A TOXIC SUBSTANCE. ALL HANDLING OF DRY MATERIALS SHOULD BE PERFORMED IN AN OPERATING FUME HOOD.

2.7.2.1 Sample Preparation--

The method of sample preparation required for XRD analysis will depend on: (1) the condition of the sample received (sample size, homogeneity, particle size distribution, and overall composition as determined by PLM); and (2) the type of XRD analysis to be performed (qualitative, quantitative, thin layer or bulk).

Bulk materials are usually received as inhomogeneous mixtures of complex composition with very wide particle size distributions. Preparation of a homogeneous, representative sample from asbestos-containing materials is particularly difficult because the fibrous nature of the asbestos minerals inhibits mechanical mixing and stirring, and because milling procedures may cause adverse lattice alterations.

A discussion of specific matrix reduction procedures is given below. Complete methods of sample preparation are detailed in Sections 2.7.2.2 and 2.7.2.3. Note: All samples should be examined microscopically before and after each matrix reduction step to monitor changes in sample particle size, composition, and crystallinity, and to ensure sample representativeness and homogeneity for analysis.

2.7.2.1.1 Milling--Mechanical milling of asbestos materials has been shown to decrease fiber crystallinity, with a resultant decrease in diffraction intensity of the specimen; the degree of lattice alteration is related to the duration and type of milling process.¹⁹⁻²² Therefore, all milling times should be kept to a minimum.

For qualitative analysis, particle size is not usually of critical importance and initial characterization of the material with a minimum of matrix reduction is often desirable to document the composition of the sample as

received. Bulk samples of very large particle size ($>2-3$ mm) should be comminuted to ~ 100 μm . A mortar and pestle can sometimes be used in size reduction of soft or loosely bound materials though this may cause matting of some samples. Such samples may be reduced by cutting with a razor blade in a mortar, or by grinding in a suitable mill (e.g., a microhammer mill or equivalent). When using a mortar for grinding or cutting, the sample should be moistened with ethanol, or some other suitable wetting agent, to minimize exposures.

For accurate, reproducible quantitative analysis, the particle size of both sample and standard materials should be reduced to ~ 10 μm (see Section 2.3.3). Dry ball milling at liquid nitrogen temperatures (e.g., Spex Freezer Mill, or equivalent) for a maximum time of 10 min is recommended to obtain satisfactory particle size distributions while protecting the integrity of the crystal lattice.⁵ Bulk samples of very large particle size may require grinding in two stages for full matrix reduction to <10 μm .^{8 16}

Final particle size distributions should always be verified by optical microscopy or another suitable method.

2.7.2.1.2 Low temperature ashing--For materials shown by PLM to contain large amounts of gypsum, cellulosic, or other organic materials; it may be desirable to ash the samples prior to analysis to reduce background radiation or matrix interference. Since chrysotile undergoes dehydroxylation at temperatures between 550°C and 650°C , with subsequent transformation to forsterite,^{23 24} ashing temperatures should be kept below 500°C . Use of a low temperature asher is recommended. In all cases, calibration of the oven is essential to ensure that a maximum ashing temperature of 500°C is not exceeded.

2.7.2.1.3 Acid leaching--Because of the interference caused by gypsum and some carbonates in the detection of asbestiform minerals by XRD (see Section 2.3.1), it may be necessary to remove these interferents by a simple acid leaching procedure prior to analysis (see Section 1.7.2.2).

2.7.2.2 Qualitative Analysis--

2.7.2.2.1 Initial screening of bulk material--Qualitative analysis should be performed on a representative, homogeneous portion of the sample with a minimum of sample treatment.

1. Grind and mix the sample with a mortar and pestle (or equivalent method, see Section 2.7.2.1.1.) to a final particle size sufficiently small ($\sim 100\ \mu\text{m}$) to allow adequate packing into the sample holder.
2. Pack the sample into a standard bulk sample holder. Care should be taken to ensure that a representative portion of the milled sample is selected for analysis. Particular care should be taken to avoid possible size segregation of the sample. (Note: Use of a back-packing method²⁵ of bulk sample preparation may reduce preferred orientation effects.)
3. Mount the sample on the diffractometer and scan over the diagnostic peak regions for the serpentine ($\sim 7.4\ \text{\AA}$) and amphibole ($8.2\text{--}8.5\ \text{\AA}$) minerals (see Table 2-2). The X-ray diffraction equipment should be optimized for intensity. A slow scanning speed of $1^\circ\ 2\theta/\text{min}$ is recommended for adequate resolution. Use of a sample spinner is recommended.
4. Submit all samples that exhibit diffraction peaks in the diagnostic regions for asbestiform minerals to a full qualitative XRD scan ($5^\circ\text{--}60^\circ\ 2\theta$; $1^\circ\ 2\theta/\text{min}$) to verify initial peak assignments and to identify potential matrix interferences when subsequent quantitative analysis is to be performed.
5. Compare the sample XRD pattern with standard reference powder diffraction patterns (i.e., JCPDS powder diffraction data³ or those of other well-characterized reference materials). Principal lattice spacings of asbestiform minerals are given in Table 2-2; common constituents of bulk insulation and wall materials are listed in Table 2-3.

2.7.2.2.2 Detection of minor or trace constituents--Routine screening of bulk materials by XRD may fail to detect small concentrations (<5 percent) of asbestos. The limits of detection will, in general, be improved if matrix absorption effects are minimized, and if the sample particle size is reduced to the optimal 1 to $10\ \mu\text{m}$ range, provided that the crystal lattice is not degraded in the milling process. Therefore, in those instances where confirmation of the presence of an asbestiform mineral at very low levels is required,

or where a negative result from initial screening of the bulk material by XRD (see Section 2.7.2.2.1) is in conflict with previous PLM results, it may be desirable to prepare the sample as described for quantitative analysis (see Section 2.7.2.3) and step-scan over appropriate 2θ ranges of selected diagnostic peaks (Table 2-2). Accurate transfer of the sample to the silver membrane filter is not necessary unless subsequent quantitative analysis is to be performed.

2.7.2.3 Quantitative Analysis--

The proposed method for quantitation of asbestos in bulk samples is a modification of the NIOSH-recommended thin-layer method for chrysotile in air.⁵ A thick-layer or bulk method involving pelletizing the sample may be used for semiquantitative analysis;^{7 8} however, this method requires the addition of an internal standard, use of a specially fabricated sample press, and relatively large amounts of standard reference materials. Additional research is required to evaluate the comparability of thin- and thick-layer methods for quantitative asbestos analysis.

For quantitative analysis by thin-layer methods, the following procedure is recommended:

1. Mill and size all or a substantial representative portion of the sample as outlined in Section 2.7.2.1.1.
2. Dry at 100° C for 2 hr; cool in a desiccator.
3. Weigh accurately to the nearest 0.01 mg.
4. Samples shown by PLM to contain large amounts of cellulosic or other organic materials, gypsum, or carbonates, should be submitted to appropriate matrix reduction procedures described in Sections 2.7.2.1.2 and 2.7.2.1.3. After ashing and/or acid treatment, repeat the drying and weighing procedures described above, and determine the percent weight loss, L.
5. Quantitatively transfer an accurately weighed amount (50-100 mg) of the sample to a 1-L volumetric flask with approximately 200 mL isopropanol to which 3 to 4 drops of surfactant have been added.
6. Ultrasonicate for 10 min at a power density of approximately 0.1 W/mL, to disperse the sample material.
7. Dilute to volume with isopropanol.

8. Place flask on a magnetic stirring plate. Stir.
9. Place a silver membrane filter on the filtration apparatus, apply a vacuum, and attach the reservoir. Release the vacuum and add several milliliters of isopropanol to the reservoir. Vigorously hand shake the asbestos suspension and immediately withdraw an aliquot from the center of the suspension so that total sample weight, W_T , on the filter will be approximately 1 mg. Do not adjust the volume in the pipet by expelling part of the suspension; if more than the desired aliquot is withdrawn, discard the aliquot and resume the procedure with a clean pipet. Transfer the aliquot to the reservoir. Filter rapidly under vacuum. Do not wash the reservoir walls. Leave the filter apparatus under vacuum until dry. Remove the reservoir, release the vacuum, and remove the filter with forceps. (Note: Water-soluble matrix interferences such as gypsum may be removed at this time by careful washing of the filtrate with distilled water. Extreme care should be taken not to disturb the sample.)
10. Attach the filter to a flat holder with a suitable adhesive and place on the diffractometer. Use of a sample spinner is recommended.
11. For each asbestos mineral to be quantitated select a reflection (or reflections) that has been shown to be free from interferences by prior PLM or qualitative XRD analysis and that can be used unambiguously as an index of the amount of material present in the sample (see Table 2-2).
12. Analyze the selected diagnostic reflection(s) by step scanning in increments of $0.02^\circ 2\theta$ for an appropriate fixed time and integrating the counts. (A fixed count scan may be used alternatively; however, the method chosen should be used consistently for all samples and standards.) An appropriate scanning interval should be selected for each peak, and background corrections made. For a fixed time scan, measure the background on each side of the peak for one-half the peak-scanning time. The net intensity, I_a , is the difference between the peak integrated count and the total background count.
13. Determine the net count, I_{Ag} , of the filter 2.36 \AA silver peak following the procedure in step 12. Remove the filter from the

holder, reverse it, and reattach it to the holder. Determine the net count for the unattenuated silver peak, I_{Ag}^o . Scan times may be less for measurement of silver peaks than for sample peaks; however, they should be constant throughout the analysis.

14. Normalize all raw, net intensities (to correct for instrument instabilities) by referencing them to an external standard (e.g., the 3.34 Å peak of an α -quartz reference crystal). After each unknown is scanned, determine the net count, I_r^o , of the reference specimen following the procedure in step 12. Determine the normalized intensities by dividing the peak intensities by I_r^o :

$$\hat{I}_a = \frac{I_a}{I_r^o}, \quad \hat{I}_{Ag} = \frac{I_{Ag}}{I_r^o}, \quad \text{and} \quad \hat{I}_{Ag}^o = \frac{I_{Ag}^o}{I_r^o}$$

2.8 CALIBRATION

2.8.1 Preparation of Calibration Standards

1. Mill and size standard asbestos materials according to the procedure outlined in Section 2.7.2.1.1. Equivalent, standardized matrix reduction and sizing techniques should be used for both standard and sample materials.
2. Dry at 100° C for 2 hr; cool in a desiccator.
3. Prepare two suspensions of each standard in isopropanol by weighing approximately 10 and 50 mg of the dry material to the nearest 0.01 mg. Quantitatively transfer each to a 1-L volumetric flask with approximately 200 mL isopropanol to which a few drops of surfactant have been added.
4. Ultrasonicate for 10 min at a power density of approximately 0.1 W/mL, to disperse the asbestos material.
5. Dilute to volume with isopropanol.
6. Place the flask on a magnetic stirring plate. Stir.
7. Prepare, in triplicate, a series of at least five standard filters to cover the desired analytical range, using appropriate aliquots of the 10 and 50 mg/L suspensions and the following procedure.

Mount a silver membrane filter on the filtration apparatus.

Place a few milliliters of isopropanol in the reservoir. Vigorously hand shake the asbestos suspension and immediately withdraw an aliquot from the center of the suspension. Do not adjust the volume in the pipet by expelling part of the suspension; if more than the desired aliquot is withdrawn, discard the aliquot and resume the procedure with a clean pipet. Transfer the aliquot to the reservoir. Keep the tip of the pipet near the surface of the isopropanol. Filter rapidly under vacuum. Do not wash the sides of the reservoir. Leave the vacuum on for a time sufficient to dry the filter. Release the vacuum and remove the filter with forceps.

2.8.2 Analysis of Calibration Standards

1. Mount each filter on a flat holder. Perform step scans on selected diagnostic reflections of the standards and reference specimen using the procedure outlined in Section 2.7.2.3, step 12, and the same conditions as those used for the samples.
2. Determine the normalized intensity for each peak measured, \hat{I}_{std}^o , as outlined in Section 2.7.2.3, step 14.

2.9 CALCULATIONS

For each asbestos reference material, calculate the exact weight deposited on each standard filter from the concentrations of the standard suspensions and aliquot volumes. Record the weight, w , of each standard. Prepare a calibration curve by regressing \hat{I}_{std}^o on w . Poor reproducibility (± 15 percent RSD) at any given level indicates problems in the sample preparation technique, and a need for new standards. The data should fit a straight line equation.

Determine the slope, m , of the calibration curve in counts/microgram. The intercept, b , of the line with the \hat{I}_{std}^o axis should be approximately zero. A large negative intercept indicates an error in determining the background. This may arise from incorrectly measuring the baseline or from interference by another phase at the angle of background measurement. A large positive intercept indicates an error in determining the baseline or that an impurity is included in the measured peak.

Using the normalized intensity, \hat{I}_{Ag} , for the attenuated silver peak of a sample, and the corresponding normalized intensity from the unattenuated

silver peak, \hat{I}_{Ag}^o , of the sample filter, calculate the transmittance, T , for each sample as follows:^{26 27}

$$T = \frac{\hat{I}_{Ag}}{I_{Ag}^o}$$

Determine the correction factor, $f(T)$, for each sample according to the formula:

$$f(T) = \frac{-R (\ln T)}{1 - T^R}$$

where

$$R = \frac{\sin \theta_{Ag}}{\sin \theta_a}$$

θ_{Ag} = angular position of the measured silver peak (from Bragg's Law), and

θ_a = angular position of the diagnostic asbestos peak.

Calculate the weight, W_a , in micrograms, of the asbestos material analyzed for in each sample, using the appropriate calibration data and absorption corrections:

$$W_a = \frac{\hat{I}_a f(t) - b}{m}$$

Calculate the percent composition, P_a , of each asbestos mineral analyzed for in the parent material, from the total sample weight, W_T , on the filter:

$$P_a = \frac{W_a (1 - 0.01L)}{W_T} \times 100$$

where

P_a = percent asbestos mineral in parent material;

W_a = mass of asbestos mineral on filter, in μg ;

W_T = total sample weight on filter, in μg ;

L = percent weight loss of parent material on ashing and/or acid treatment (see Section 2.7.2.3).

2.10 REFERENCES

1. H. P. Klug and L. E. Alexander, X-ray Diffraction Procedures for Polycrystalline and Amorphous Materials, 2nd ed., New York: John Wiley and Sons, 1979.
2. L. V. Azaroff and M. J. Buerger, The Powder Method of X-ray Crystallography, New York: McGraw-Hill, 1958.
3. JCPDS-International Center for Diffraction Data Powder Diffraction File, U.S. Department of Commerce, National Bureau of Standards, and Joint Committee on Powder Diffraction Studies, Swarthmore, PA.
4. W. J. Campbell, C. W. Huggins, and A. G. Wylie, Chemical and Physical Characterization of Amosite, Chrysotile, Crocidolite, and Nonfibrous Tremolite for National Institute of Environmental Health Sciences Oral Ingestion Studies, U.S. Bureau of Mines Report of Investigation RI8452, 1980.
5. B. A. Lange and J. C. Haartz, Determination of microgram quantities of asbestos by X-ray diffraction: Chrysotile in thin dust layers of matrix material, Anal. Chem., 51(4):520-525, 1979.
6. NIOSH Manual of Analytical Methods, Volume 5, U.S. Dept. HEW, August 1979, pp. 309-1 to 309-9.
7. H. Dunn and J. H. Stewart, Jr., Quantitative determination of chrysotile in building materials, The Microscope, 29(1), 1981.
8. M. Taylor, Methods for the quantitative determination of asbestos and quartz in bulk samples using X-ray diffraction, The Analyst, 103(1231): 1009-1020, 1978.
9. L. Birks, M. Fatemi, J. V. Gilfrich, and E. T. Johnson, Quantitative Analysis of Airborne Asbestos by X-ray Diffraction, Naval Research Laboratory Report 7879, Naval Research Laboratory, Washington, D.C., 1975.
10. U.S. Environmental Protection Agency, Asbestos-Containing Materials in School Buildings: A Guidance Document, Parts 1 and 2, EPA/OTS No. C00090, March 1979.
11. J. B. Krause and W. H. Ashton, Misidentification of asbestos in talc, pp. 339-353, in: Proceedings of Workshop on Asbestos: Definitions and Measurement Methods (NBS Special Publication 505), C. C. Gravatt, P. D. LaFleur, and K. F. Heinrich (eds.), Washington, D.C.: National Measurement Laboratory, National Bureau of Standards, 1977 (issued 1978).
12. H. D. Stanley, The detection and identification of asbestos and asbestiform minerals in talc, pp. 325-337, in Proceedings of Workshop on Asbestos: Definitions and Measurement Methods (NBS Special Publication 506), C. C. Gravatt, P. D. LaFleur, and K. F. Heinrich (eds.), Washington, D.C.

National Measurement Laboratory, National Bureau of Standards, 1977 (issued 1978).

13. A. L. Rickards, Estimation of trace amounts of chrysotile asbestos by X-ray diffraction, Anal. Chem., 44(11):1872-3, 1972.
14. P. M. Cook, P. L. Smith, and D. G. Wilson, Amphibole fiber concentration and determination for a series of community air samples: use of X-ray diffraction to supplement electron microscope analysis, in: Electron Microscopy and X-Ray Applications to Environmental and Occupation Health Analysis, P. A. Russell and A. E. Hutchings (eds.), Ann Arbor: Ann Arbor Science Publications, 1977.
15. A. N. Rohl and A. M. Langer, Identification and quantitation of asbestos in talc, Environ. Health Perspectives, 9:95-109, 1974.
16. J. L. Graf, P. K. Ase, and R. G. Draftz, Preparation and Characterization of Analytical Reference Minerals, DHEW (NIOSH) Publication No. 79-139, June 1979.
17. J. C. Haartz, B. A. Lange, R. G. Draftz, and R. F. Scholl, Selection and characterization of fibrous and nonfibrous amphiboles for analytical methods development, pp. 295-312, in: Proceedings of Workshop on Asbestos: Definitions and Measurement Methods (NBS Special Publication 506), C. C. Gravatt, P. D. LaFleur, and K. F. Heinrich (eds.), Washington, D.C.: National Measurement Laboratory, National Bureau of Standards, 1977 (issued 1978).
18. Personal communication, A. M. Langer, Environmental Sciences Laboratory, Mount Sinai School of Medicine of the City University of New York, New York, New York.
19. A. M. Langer, M. S. Wolff, A. N. Rohl, and I. J. Selikoff, Variation of properties of chrysotile asbestos subjected to milling, J. Toxicol. and Environ. Health, 4:173-188, 1978.
20. A. M. Langer, A. D. Mackler, and F. D. Pooley, Electron microscopical investigation of asbestos fibers, Environ. Health Perspect., 9:63-80, 1974.
21. E. Occella and G. Maddalon, X-ray diffraction characteristics of some types of asbestos in relation to different techniques of comminution, Med. Lavoro, 54(10):628-636, 1963.
22. K. R. Spurny, W. Stöber, H. Opiela, and G. Weiss, On the problem of milling and ultrasonic treatment of asbestos and glass fibers in biological and analytical applications, Am. Ind. Hyg. Assoc. J., 41:198-203, 1980.
23. L. G. Berry and B. Mason, Mineralogy, San Francisco: W. H. Greeman & Co., 1959.

24. J. P. Schelz, The detection of chrysotile asbestos at low levels in talc by differential thermal analysis, Thermochimica Acta, 8:197-204, 1974.
25. Reference 1, pp. 372-374.
26. J. Leroux, Staub-Reinhalt Luft, 29:26 (English), 1969.
27. J. A. Leroux, B. C. Davey, and A. Paillard, Am. Ind. Hyg. Assoc. J., 34:409, 1973.

